


# NORLAND PRODUCTS INCORPORATED


2540 Route 130, Building 100, Cranbury, NJ  
Mailing Address • P.O. Box 637 • Cranbury, NJ 08512  
Tel • 609-395-196 Fax • 609-395-9006

## Shipping Kit Instructions

1. Place the shipping block on the right side of the interferometer with the two screws and external washers as shown in figure 1. It may be necessary to adjust the X & Y knobs to line up the holes for the shipping block.

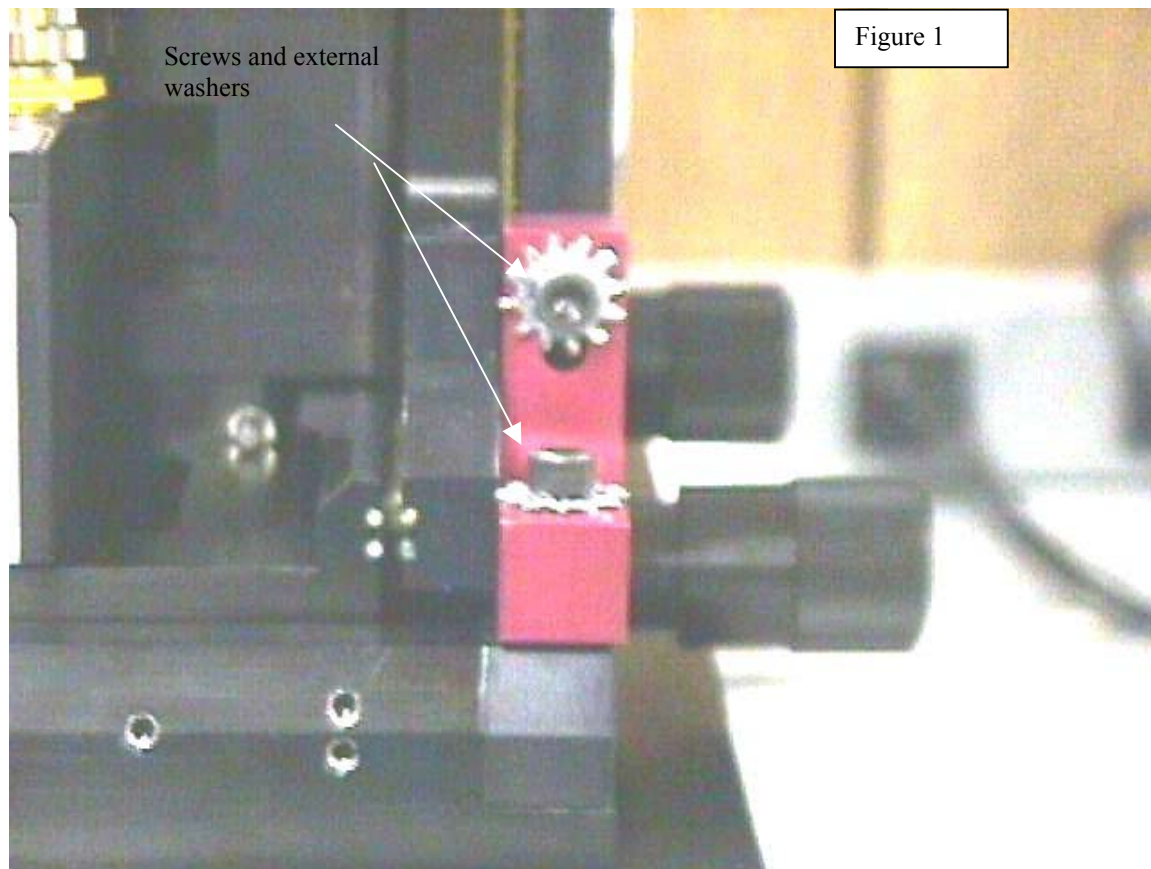


2. Place the shipping screw into the Tilt Stage as shown in figure 2.

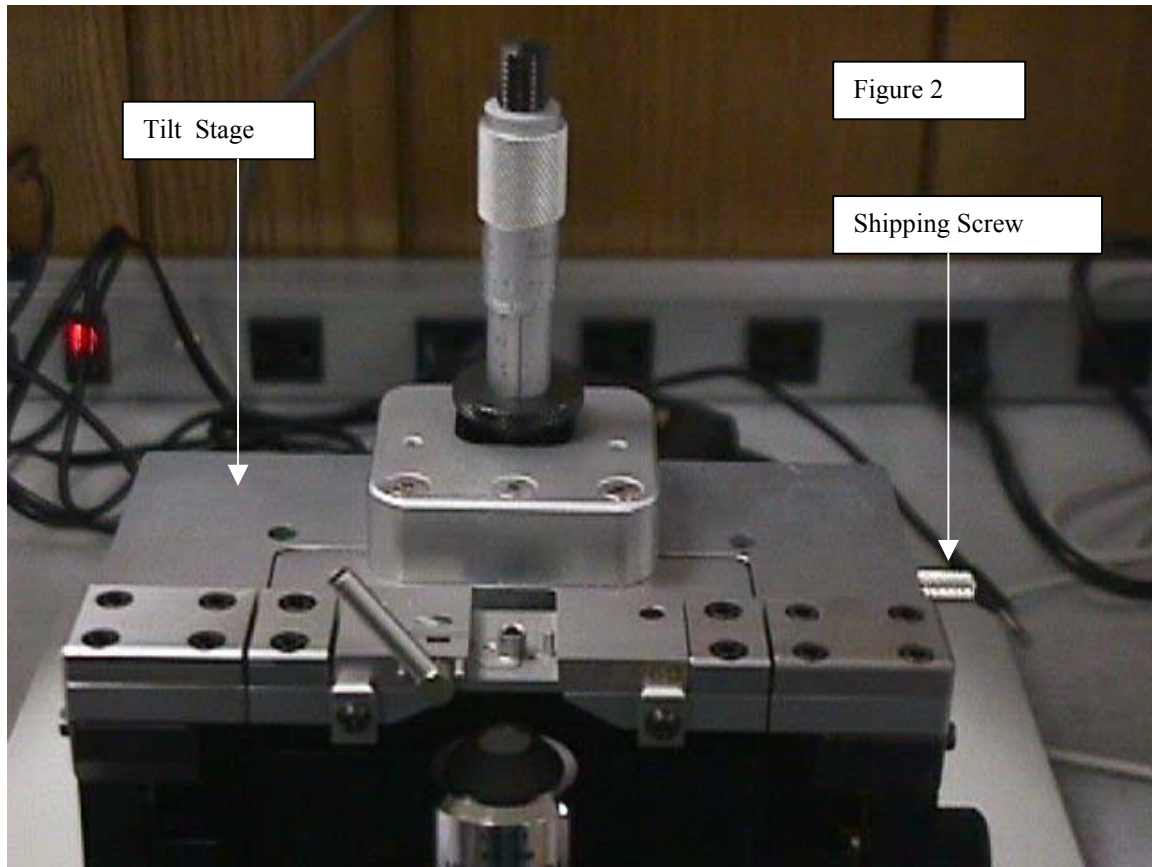

3. Remove the Tilt Stage from the interferometer by removing the four screws and place it into the accessories box as shown in figure 3. Add foam to protect the Tilt Stage.




- Next place the Objective Insert over the 10X lens as shown in figure 4.

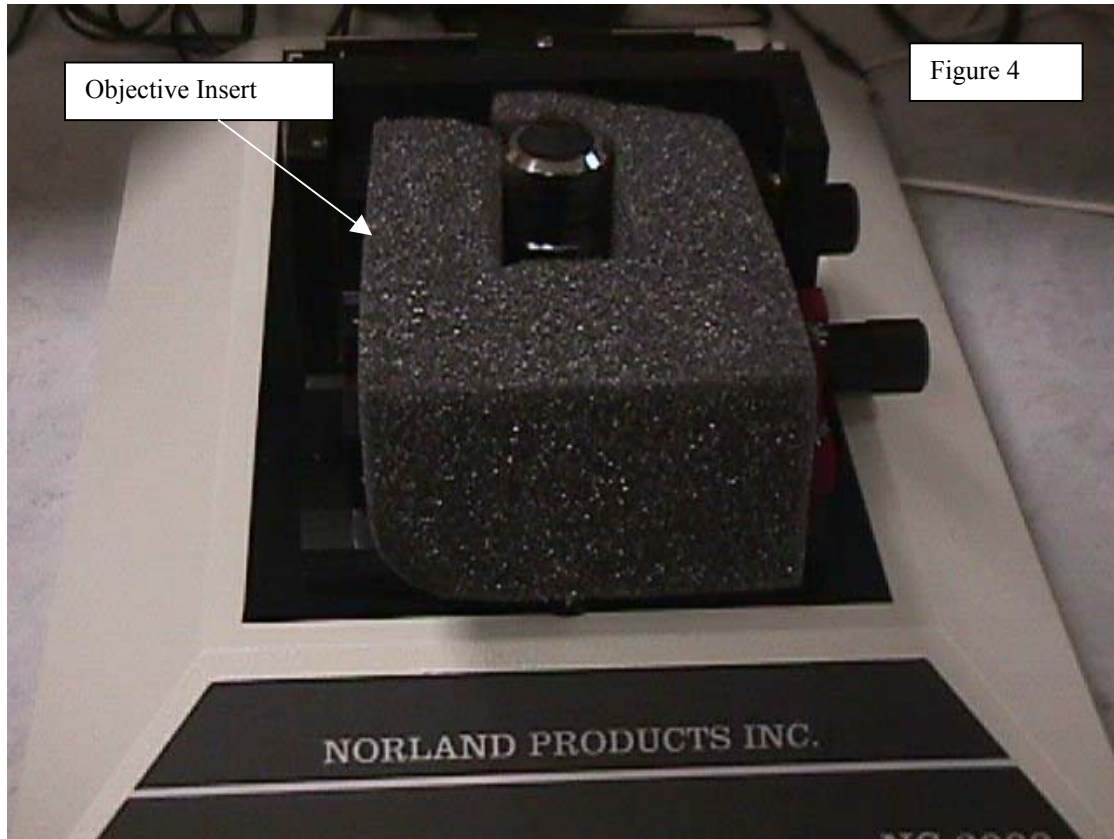

- Next place the interferometer into the plastic bag as shown in figure 5.



6. Prepare the shipping box and place the microscope insert into the box as shown in figure 6.

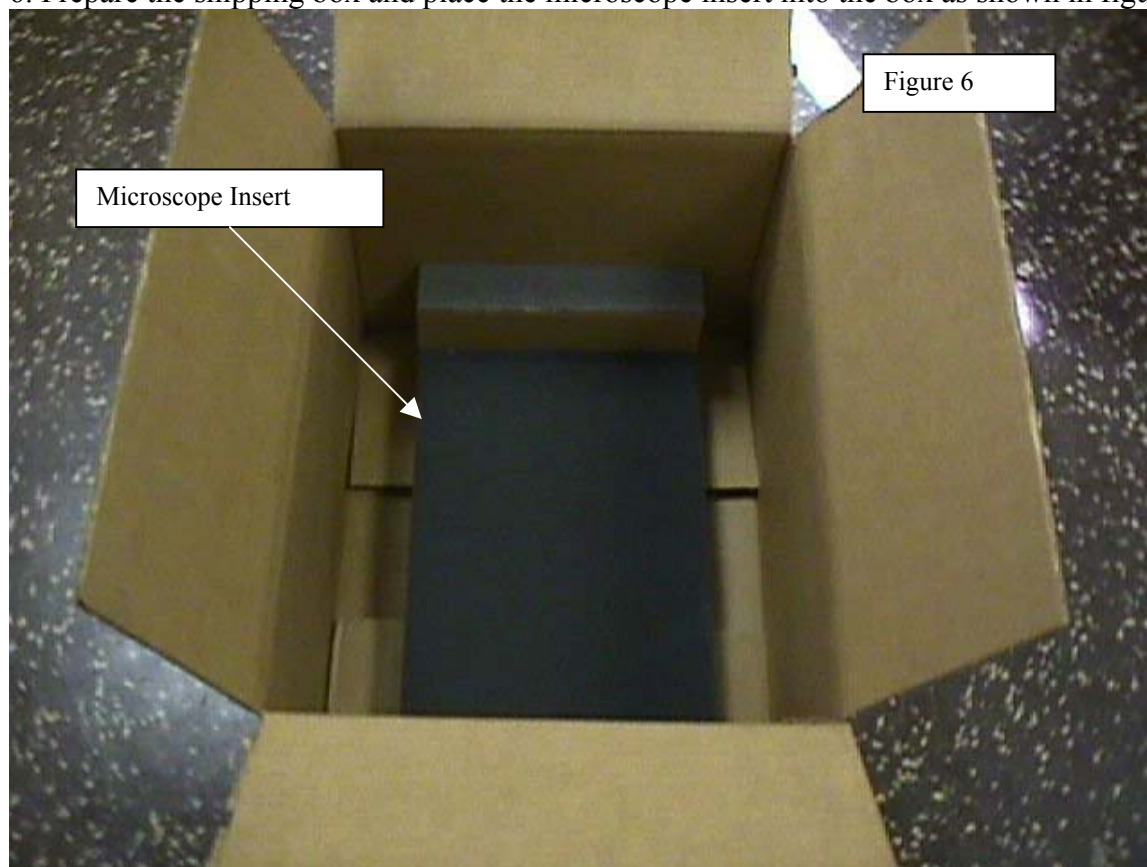


6.

7. Proceed to place the interferometer into the shipping box as shown in figure 7.




8. Place the right and left inserts into the shipping box as shown in figure 8.

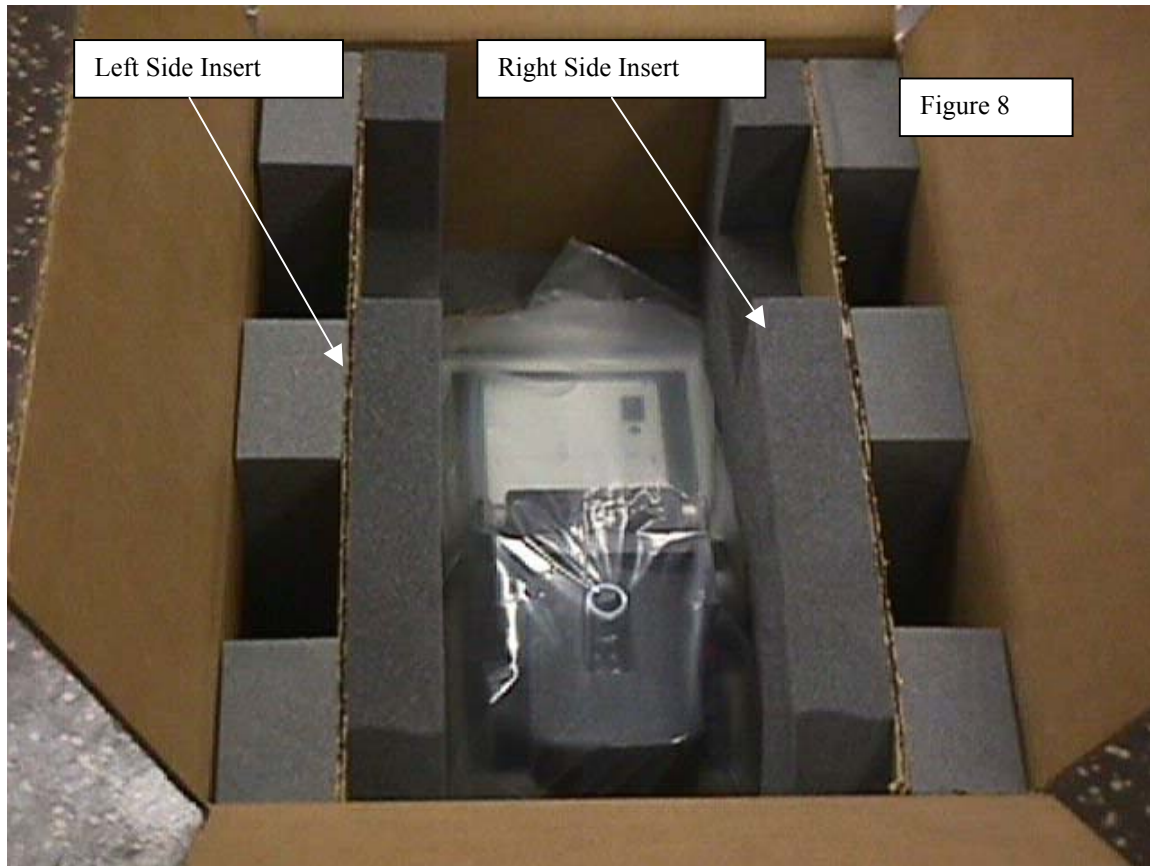

9. Place the accessories box into the shipping box as shown in figure 9.



10. Place the microscope spacer over the 10x lens as shown in figure 10.

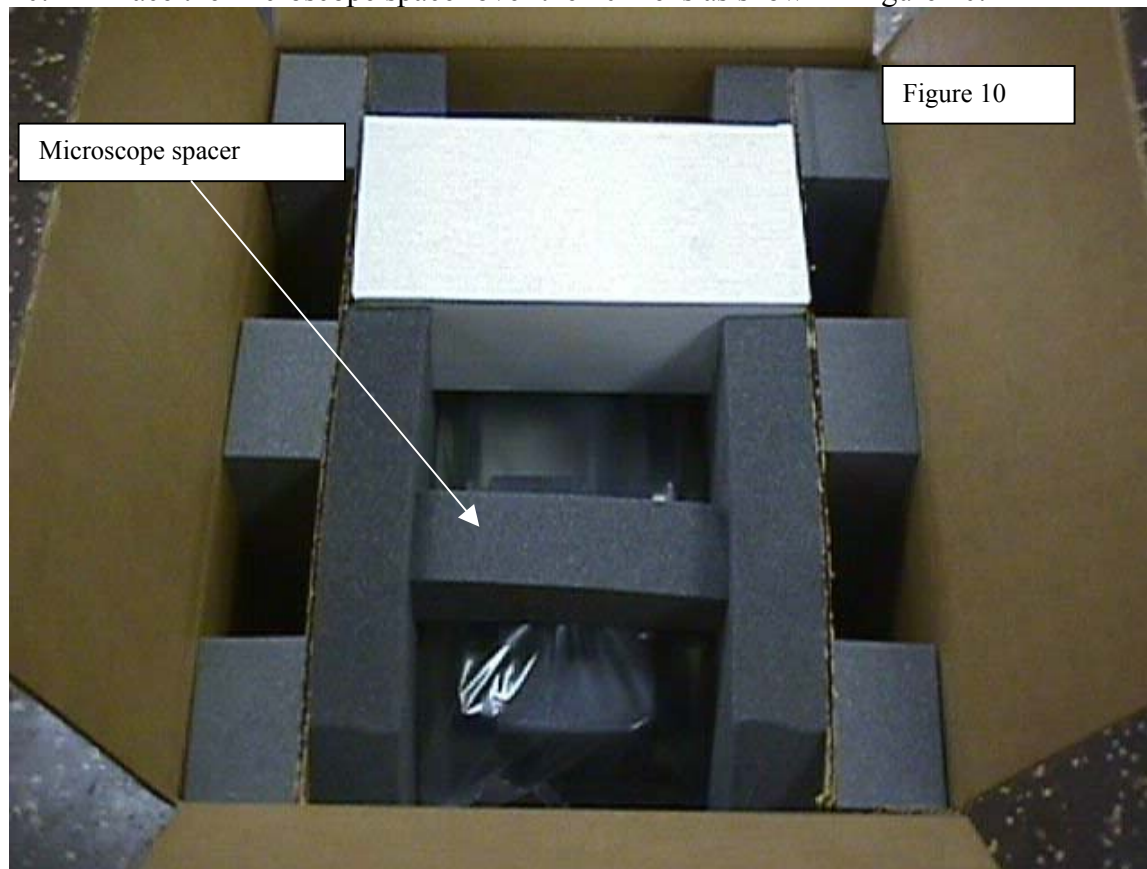

11. Close the box and seal it with strong shipping tape as shown in figure 11.

